

# STANDARD RESPONSE LEVER

ENV-110M, ENV-110RM, ENV-116M, ENV-610

## WIRING GUIDE


### STANDARD RESPONSE LEVER

When the lever is pressed, the micro-switch is closed. The closing of the switch causes the input bit to be pulled (connected) to ground signaling a response from the lever at the computer.

### WIRING

The input cable connects to any available Input on a Standard MED Connection Panel.

Molex Pin #	Function	Wire Color
1	28V Ground	Black
2	Lever Response	Red or White
3	Not Used	

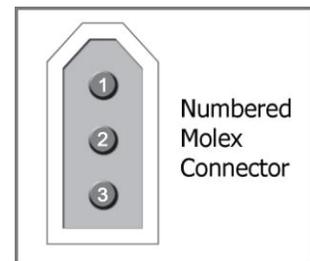

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at [www.med-associates.com](http://www.med-associates.com) for contact information.

For technical questions, email [support@med-associates.com](mailto:support@med-associates.com).

DOC-045

Rev. 1.1

Copyright ©2012  
All Rights Reserved

Med Associates Inc.  
P.O. Box 319  
St. Albans, Vermont 05478

Phone: 802.527.2343  
Fax: 802.527.5095  
[www.med-associates.com](http://www.med-associates.com)