

RESPONSE FEEDBACK RELAY

ENV-135M

WIRING GUIDE

DOC-061

Rev. 1.1

Copyright ©2012

All Rights Reserved

Med Associates Inc.

P.O. Box 319

St. Albans, Vermont 05478

Phone: 802.527.2343

Fax: 802.527.5095


www.med-associates.com

CHAPTER 1 | WIRING

OPERATE CABLE

The ENV-135M Cable should be connected to any available Output on a standard MED Connection Panel. The relay will click when the corresponding Output bit is turned “ON”. Also, by physically connecting the white wire to ground the relay will click.

Molex Pin #	Function	Wire Color
1	Not Used	
2	Operate	White
3	+28 Volts	Red

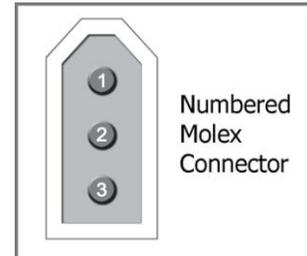

APPENDIX A | CONTACT INFORMATION

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.