

TRIPLE STIMULUS LIGHT

ENV-222 & ENV-322

HARDWARE MANUAL

DOC-067

Rev. 1.2

Copyright ©2013
All Rights Reserved

Med Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478


Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

CHAPTER 1 | WIRING

The Triple Stimulus Light includes cable part number SG-223. This cable terminates with three mini Molex connectors. Each of these connectors should be connected to any available Output ports on a Standard MED Connection Panel. Each LED will illuminate when the corresponding output bit is turned “ON”.

NOTE: The connector that controls the yellow LED (indicated in the diagram below) supplies power to each of the three LEDs, and therefore must be connected even if the yellow LED is not being used.

If a connection panel is not being used, connecting the red wire to +28 Volts, and the black, white, or green wire to 28 Volt ground will cause the corresponding LED to illuminate.

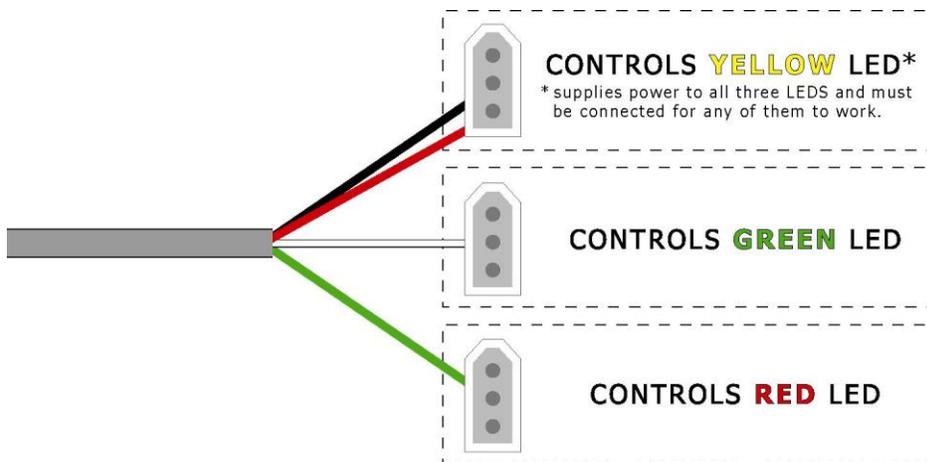

APPENDIX A | CONTACT INFORMATION

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.