

LIGHTS

ENV-315M HOUSE LIGHT

ENV-321M STIMULUS LIGHT

WIRING GUIDE

DOC-081

Rev. 1.2

Copyright ©2012

All Rights Reserved

Med Associates Inc.

P.O. Box 319

St. Albans, Vermont 05478

Phone: 802.527.2343

Fax: 802.527.5095


www.med-associates.com

CHAPTER 1 | WIRING

Output Cable

Connect the light's output cable to any available Output on a Standard MED Connection Panel. The light will turn on when a signal is sent to the output.

Molex Pin #	Function	Wire Color
1	Not Used	
2	Light Activate	White
3	+28 Volts	Red

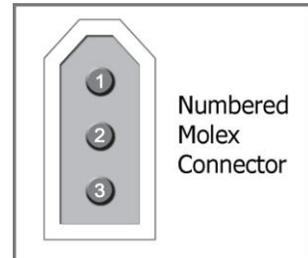

APPENDIX A | CONTACT INFORMATION

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.