

BARNES MAZE

ASSEMBLY INSTRUCTIONS



ENV-562 Users Manual
DOC-098
Rev. 1.2

Copyright © 2013
All Rights Reserved

MED Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478
www.med-associates.com

Hardware Guide

The Barnes Maze includes the following hardware:

- Base Support
- Center Column
- Maze Top
- Goal Box
- Eight Retaining Pins

Assembly Instructions

1. Place the base support on the floor with the alignment pins up, as shown in Figure 1.

Figure 1 – Properly Oriented Base Support



2. Orient the center column so that the alignment pins are up. Connect the center column to the base as shown in Figure 2.

Figure 2 – Base Support and Center Column Assembly



3. Secure the center column to the base support by inserting a retaining pin into each of the four holes around the bottom of the center column, as shown in Figure 3.

Figure 3 – Secure Center Column with Retaining Pins



4. Connect the maze top to the center column as shown in Figure 4.

Figure 4 – Maze Top Placed on the Center Column



5. Secure the maze top to the center column by inserting a retaining pin into each of the four holes around the top of the center column, as shown in Figure 5.

Figure 5 – Maze Top Secured to the Center Column



6. Install the goal box by sliding it into the track under the desired hole.

Figure 6 – Installed Goal Box

