

PIGEON KEY W/MULTI-COLOR DISPLAY

ENV-124AM

USER'S MANUAL

DOC-134

Rev. 1.0

Copyright ©2014
All Rights Reserved

Med Associates, Inc.
P.O. Box 319
St. Albans, Vermont 05478

Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

LIGHT CABLE

The Light Cable terminates with 12 Molex connectors that should be connected to any available Outputs on a Standard MED Connection Panel. Each connector controls a different color light, as shown in the table below. Note that Connector 1 must be connected in order for any other light to function.

Connector #	Molex Pin #				
	1	Wire Color	Display Color	Wire Color	Function
1	Not Used	White	White	Red	+ 28 Volts
2	Not Used	Black	Red 1	Not Used	
3	Not Used	Green	Red 2	Not Used	
4	Not Used	Blue	Red 3	Not Used	
5	Not Used	Yellow	Amber	Not Used	
6	Not Used	Orange	Yellow	Not Used	
7	Not Used	Brown	Green 1	Not Used	
8	Not Used	Gray	Green 2	Not Used	
9	Not Used	Pink	Blue 1	Not Used	
10	Not Used	Violet	Blue 2	Not Used	
11	Not Used	Tan	Blue 3	Not Used	
12	Not Used	Red/Black	Blue 4	Not Used	

RESPONSE CABLE

Molex Pin #	Wire Color	Function
1	Black	28 Volt Ground
2	White	Key Response
3	Not Used	

