

RESPONSE WHEEL


ENV-113M

USER'S GUIDE

CONNECTIONS

The Response Wheel cable should be connected to any available Input on a Standard Med connection panel. A switch closure is produced every 90 degrees of rotation.

Molex Pin #	Function	Wire Color
1	28V Ground	Black
2	Input Bit	White
3	Not Used	

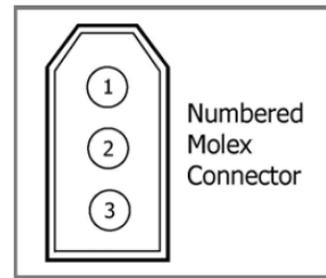

INSTALLATION

1. Remove the two screws connecting the main wheel panel to the modular panel.
2. Slide the modular panel into the channels of one bay in the chamber.
3. Attach the main wheel to the modular panel using the two screws removed above.

DOC-136

Rev. 1.0

Copyright ©2014
All Rights Reserved

Med Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478

Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com