

MED-PC[®] IV USB INTERFACE

INSTALLATION MANUAL FOR MED-PC[®] WORKSHOP

DOC-138
Rev. 1.0

Copyright © 2007
All Rights Reserved

MED Associates, Inc.
P.O. Box 319
St. Albans, Vermont 05478
www.med-associates.com

INTRODUCTION


Thank you for attending the MED-PC® Training Session. This course will use MED-PC® software in conjunction with our convenient DIG-703 USB Interface. This manual describes the installation procedure that will be used to install both the MED-PC® IV software and the DIG-703 USB Interface drivers on the laptop computer.

NOTE: Please do not connect the DIG-703 Interface to the computer until after the software has been installed.

SOFTWARE INSTALLATION

Begin by inserting the MED-PC® IV USB Interface Installation CD into the CD-ROM drive. The screen shown in Figure 1 will appear.

Figure 1 – Main Screen

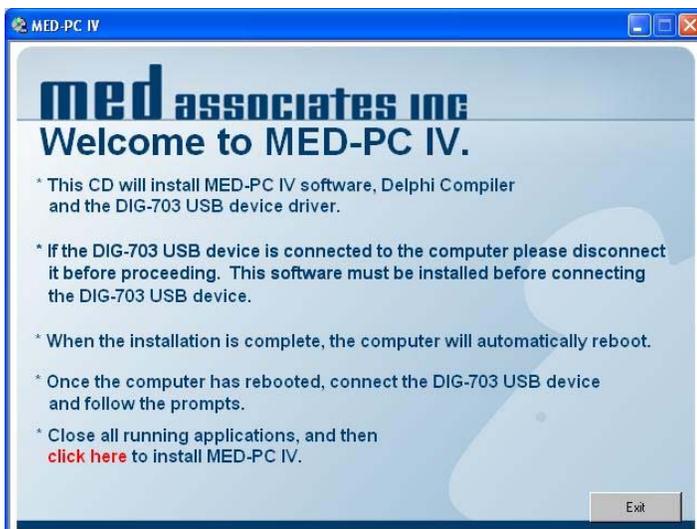

As stated on the screen above, verify that the DIG-703A USB Interface is not connected to the computer, and then close all running applications. Next, click on “click here” to begin installation. The MED-PC® software and all the related files that are necessary, such as sample MedState® notation files and DIG-703 USB drivers, will automatically be installed. This procedure will take several minutes. During this time, a brief overview of the MED Associates product catalog will be displayed. When installation is completed, the computer will automatically reboot.

DIG-703A USB INTERFACE INSTALLATION

Once the computer has rebooted, use the included power cord to plug the DIG-703A USB Interface in to a standard wall outlet.

Figure 2 - DIG-703A USB Interface



Figure 3 - Plug In the DIG-703A USB Interface



Be sure that the DIG-703A USB Interface is switched off, and then use the included USB cable to connect the USB port on the back of the DIG-703A to any available USB port on the computer.

Figure 4 - Connect the DIG-703A to Computer USB Port



Now that the DIG-703A USB Interface is properly connected, it may be switched on. Once the unit is turned on the Found New Hardware Wizard, shown in Figure 5, will appear. Select **Install the software automatically (Recommended)** and then click **Next**.

Figure 5 - Driver Installation



After a few seconds, the screen shown in Figure 6 will appear. The DIG-703A USB Interface driver installation is now complete.

Figure 6 - Driver Installation Complete

