

MONKEY HAND HOLE

ENV-600-11S

WIRING GUIDE

DOC-152

Rev. 1.1

Copyright ©2013

All Rights Reserved

Med Associates Inc.

P.O. Box 319

St. Albans, Vermont 05478

Phone: 802.527.2343

Fax: 802.527.5095


www.med-associates.com

CHAPTER 1 | WIRING

STIMULUS LIGHT CABLE

The Stimulus Light Cable connects to any available Output on a Standard MED Connection Panel. Turning on the corresponding output will turn on the stimulus light.

Molex Pin #	Function	Wire Color
1		Not Used
2	Stimulus Light	White
3	+28 Volts	Red

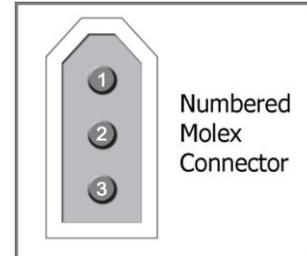

RESPONSE CABLE

The Response Cables connect to any available Inputs on a Standard MED Connection Panel.

Connector 1

Molex Pin #	Function	Wire Color
1	28V Ground	Black
2	Right Pull Switch	White
3	Not Used	

Connector 2

Molex Pin #	Function	Wire Color
1	28V Ground	Black
2	Left Pull Switch	White
3	Not Used	

APPENDIX A | CONTACT INFORMATION

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.