

MOUSE SWIVEL ARM

PHM-124M

USER'S MANUAL

DOC-160

Rev. 1.1


Copyright ©2012
All Rights Reserved

Med Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478


Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

CHAPTER 1 | ASSEMBLY

Using the ENV-307A Mouse Chamber

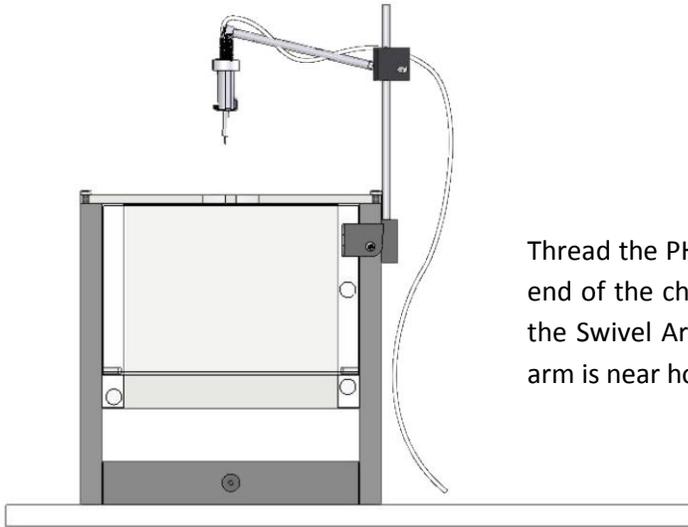


Thread the PHM-124M Swivel Arm into the block on the end of the chamber, as shown at left. Use the knob on the Swivel Arm to adjust the spring tension so that the arm is near horizontal.

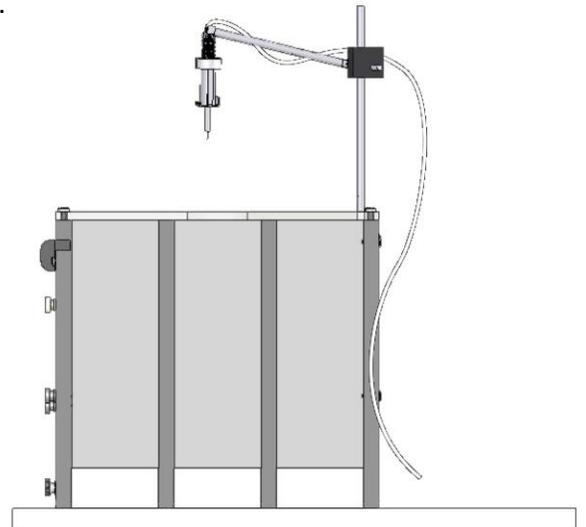

Using the ENV-307W Mouse Chamber

Thread the PHM-124M Swivel Arm into the hole on the top of the chamber, as shown at right. Use the knob on the Swivel Arm to adjust the spring tension so that the arm is near horizontal.

Inserting the Swivel

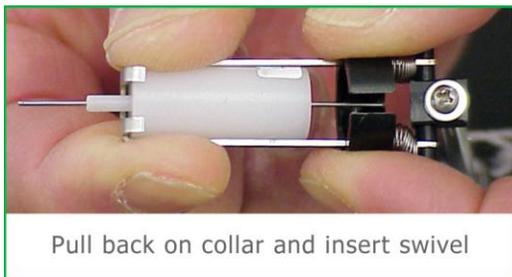

APPENDIX A | CONTACT INFORMATION

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.