

RETRACTABLE RESPONSE LEVER

ENV-112BM

WIRING GUIDE

DOC-168

Rev. 1.1

Copyright ©2013

All Rights Reserved

Med Associates Inc.

P.O. Box 319

St. Albans, Vermont 05478

Phone: 802.527.2343

Fax: 802.527.5095


www.med-associates.com

CHAPTER 1 | WIRING

OPERATE CABLE

The Operate Cable should be connected to any available Output on a standard MED Connection Panel. The +28 Volt DC motor is activated by turning the output bit “ON”, causing the lever to extend into the chamber. When the bit is turned “OFF” the lever will then retract out of the chamber.

Molex Pin #	Function	Wire Color
1	Not Used	
2	Lever Operate	White
3	+28 Volts DC	Red

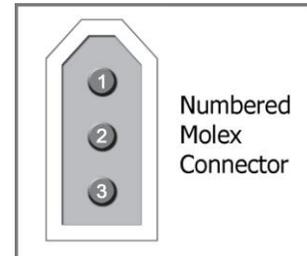

RESPONSE CABLE

The Response Cable should be connected to any available Input on a standard MED Connection Panel. When the animal depresses the lever, the signal is received via the input bit.

Molex Pin #	Function	Wire Color
1	28V Ground	Black
2	Lever Response	White
3	Not Used	

APPENDIX A | CONTACT INFORMATION

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.