


PHM-119B RAT SELF INFUSION PACKAGE

HARDWARE SHEET

This package includes one of each of the following:

MED Part Number	Description
PHM-119-1	Quick Connect™ Harness
PHM-119-2	Quick Connect™ Swivel
PHM-119-3B	12-inch Quick Connect™ Tether
PHM-119-4B	Tapered Femoral Vein Catheter

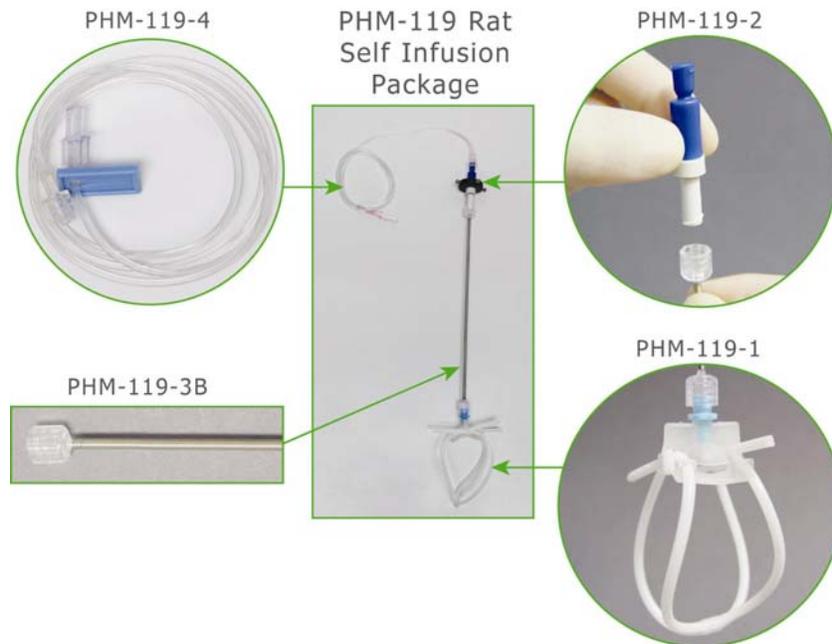

Place the Gimbal Ring from the PHM-110 Arm Assembly over the PHM-119-2 and tighten the setscrews. Thread the blue end of the PHM-119-2 onto the male end of the PHM-119-4. Next thread the white end of the PHM-119-2 onto the PHM-119-3B. Connect the PHM-119-1 to the other end of the PHM-119-3B. The completed assembly is shown above.

DOC-171
Rev. 1.0

Copyright © 2007
All Rights Reserved

MED Associates, Inc.
P.O. Box 319
St. Albans, Vermont 05478
www.med-associates.com