


ENV-130M TWELVE STIMULUS PROJECTOR WITH PECKING KEY

Pecking Key Response Cable

Connect the Response Cable to any available input on a standard MED Associates connection panel. When the key is pressed, an output will be generated.

Molex Pin #	Function	Wire Color
1	28 Volt Ground	Black
2	Key Response	Red
3	-	-

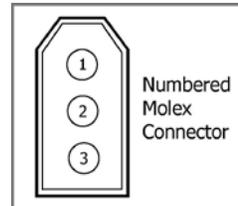

Control Cable

The control cable terminates with twelve 3-pin Molex connectors. Connector 1 (Black & Red/Green wires) as well as the connectors that correspond to the desired stimulus projections must be connected to any available output on a standard MED connection panel. Connector 1 supplies +28 Volts to the Stimulus Projector.

Connector	Stimulus Projection	Molex Pin #		
		1	2	3
1		-	Black	Red/Green
2		-	Brown	-
3		-	Red	-
4		-	Orange	-
5		-	Yellow	-
6		-	Green	-

Connector	Stimulus Projection	Molex Pin #		
		1	2	3
7		-	Blue	-
8		-	Violet	-
9		-	Gray	-
10		-	White	-
11		-	Tan	-
12		-	Pink	-

DOC-172
Rev. 1.0

Copyright © 2016
All Rights Reserved

Med Associates, Inc.
P.O. Box 319
St. Albans, Vermont 05478
www.med-associates.com