

STIMULUS LIGHT FADER CONTROL

ENV-226

USER'S MANUAL


DOC-174

Rev. 1.1

Copyright ©2014
All Rights Reserved


Med Associates, Inc.
P.O. Box 319
St. Albans, Vermont 05478

Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

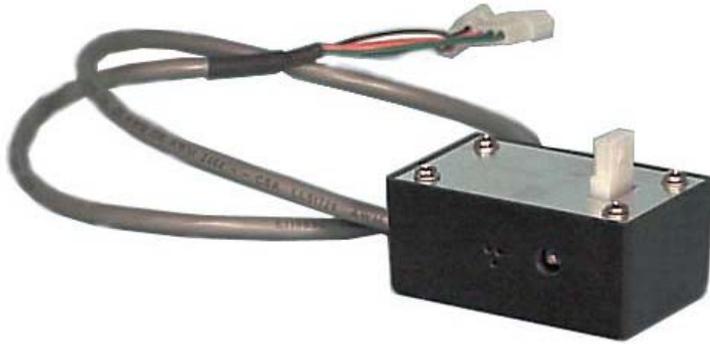

Intensity Control Cable

Molex Pin #	Connector 1		Connector 2	
	Function	Wire Color	Function	Wire Color
1	28V Ground	Black	28V Ground	Green
2	Full Intensity	White	Dimmer Activate	Blue
3	+28 Volts	Red	+28 Volts	Orange

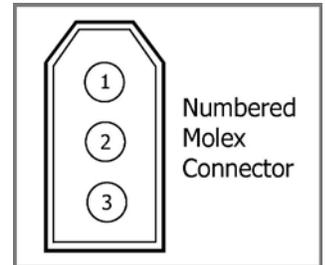

The ENV-226 Stimulus Light Fader control can be used with the stimulus light (ENV-221), the house light (ENV-215M) or stimulus/house light (ENV-227M). Connect the light cable to the 3-pin male Molex connector on the top of the ENV-226 control box.

Connectors 1 and 2 should be connected to any available outputs on a standard Med Associates connection panel. To operate the light at full intensity the output bit controlling Connector 1 must be turned “ON”. For less than full intensity the output bit controlling Connector 2 must be turned “ON”. If both output bits are turned “ON” the full intensity will override the dimmed intensity.

The intensity is adjusted using the dimmer control on the side of the black control box. The intensity is factory set to approximately half of the full intensity (15 Volts on the output connector to the light).

CONTACT INFORMATION

Please contact Med Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.