

STARTLE PLATFORM

PHM-250

LOAD CELL ASSEMBLY INSTRUCTIONS

DOC-219

Rev. 1.1

Copyright ©2013
All Rights Reserved

Med Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478

Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

CHAPTER 1 | INTRODUCTION

The PHM-250 Load Cell is shipped disassembled in order to protect the sensitive hardware. Consequently, some assembly is required.

Figure 1-1 – Load Cell Platform, unassembled




To assemble the platform, align the Load Cell housing over the base and press into place. The housing should be oriented so that the cable is on the right side. Secure the housing in place using the latches on the opposite end (see Figure 1-2).

Figure 1-2 – Align housing over the base



Place the Load Cell Platform over the housing. Orient the platform so that the groove in the bottom is aligned with the load cell, as shown in Figure 1-3.

Figure 1-3 – Place Load Cell Platform over housing.

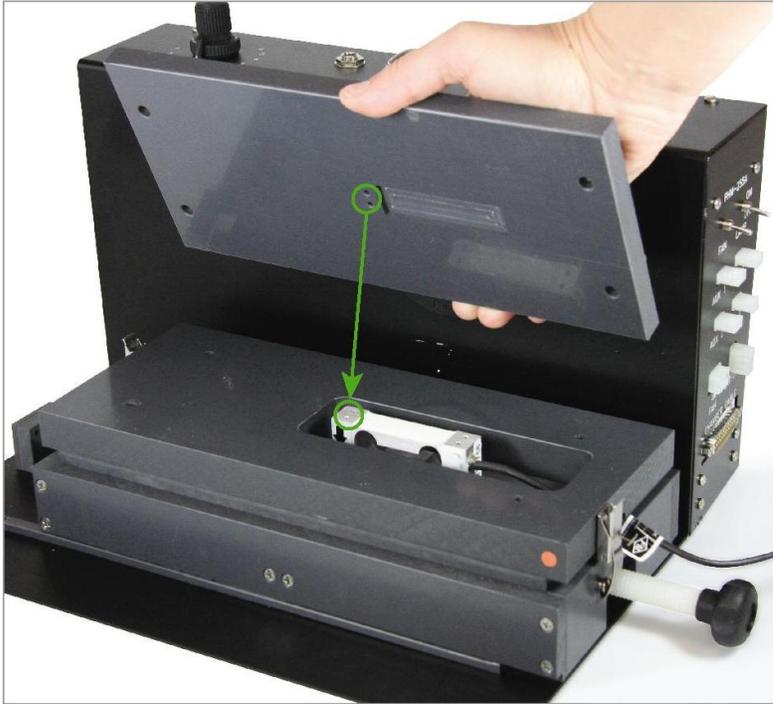

Secure the platform in place using the cap head screws (included) and a 2.5mm Allen wrench. Ensure the platform is aligned with the housing before tightening the screws.

Figure 1-4 – Secure platform with cap head screws



APPENDIX A | CONTACT INFORMATION

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.