

ENV-596M-C / ENV-596R-C

WATER MAZE POOL SUPPORT WITH CASTERS

User's Sheet

DOC-231
Rev. 1.1

Copyright ©2013
All Rights Reserved

Med Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478

Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

WATER MAZE POOL SUPPORT WITH CASTERS

The Water Maze Pool Support with Casters is equipped with levelers, allowing the user to easily relocate the Water Maze hardware.

NOTE: Levelers must be lowered prior to filling the Water Maze Pool with water. Do not attempt to move the pool while full. If the pool is equipped with a heater an/or On Demand Platform, be sure to turn off these devices before draining the pool.

Relocating the Water Maze Hardware

1. Drain the water maze pool.
2. Raise the levelers using the supplied wrench as shown in the diagram below, until the casters are touching the floor.
3. Move the empty pool and Pool Support to the desired location.
4. Lower the levelers using the supplied wrench according to the diagram below, until the casters are no longer touching the floor.
5. Observe the bubble levels to verify that the Pool Support is level.
6. Once the Pool Support wheels are no longer in contact with the floor and the Pool Support is level, the Water Maze Pool may be filled.

