

# **NIR BACKLIT MAZE**

## **ASSEMBLY MANUAL**

**DOC-243**

**Rev. 1.1**

Copyright ©2013  
All Rights Reserved

Med Associates Inc.  
P.O. Box 319  
St. Albans, Vermont 05478

Phone: 802.527.2343  
Fax: 802.527.5095  
[www.med-associates.com](http://www.med-associates.com)




## **Table of Contents**

Assembly Instructions .....	1
Disassembly Instructions .....	3
Appendix A.....	3
Contact Information.....	3

## ASSEMBLY INSTRUCTIONS


1. The first step is to identify the power runway, shown in Figure 1. This runway has a power switch on the end. There is one power runway per maze.

*Figure 1 – End of the Power Runway*

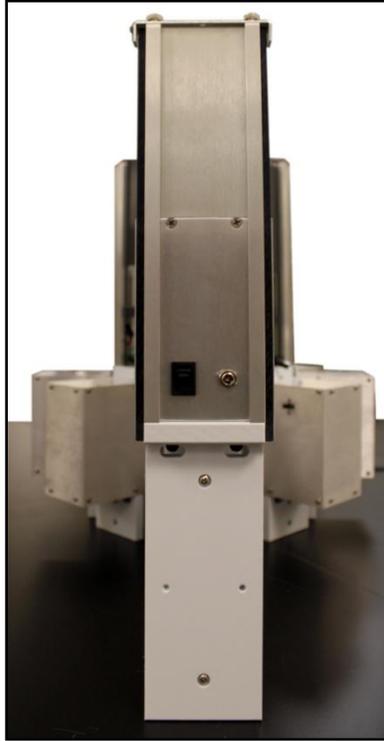


2. Next, identify the power location on the maze hub, shown in Figure 2. There is one power location per hub. This location has three wires, whereas the others have only one.

*Figure 2 - Power Location on the Hub*

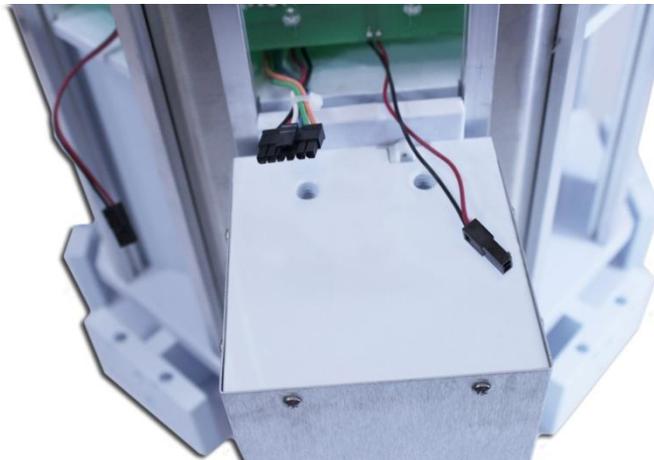


3. If not already installed on the hub, slide doors the doors into place.
4. Place the power runway close to the power location and connect the three cables, as shown in Figure 3.

*Figure 3 - Cables Connected*

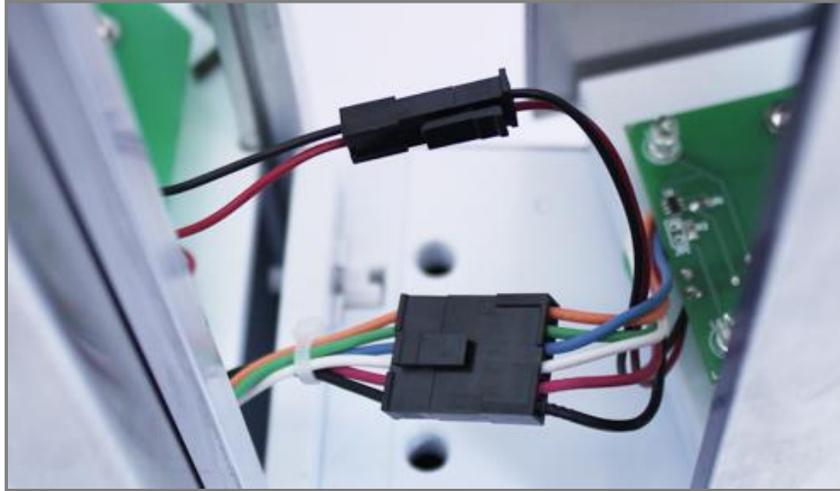

5. Tuck the connectors and excess cable underneath the circuit boards, as shown in Figure 4.

*Figure 4 - Cables under Circuit Board*

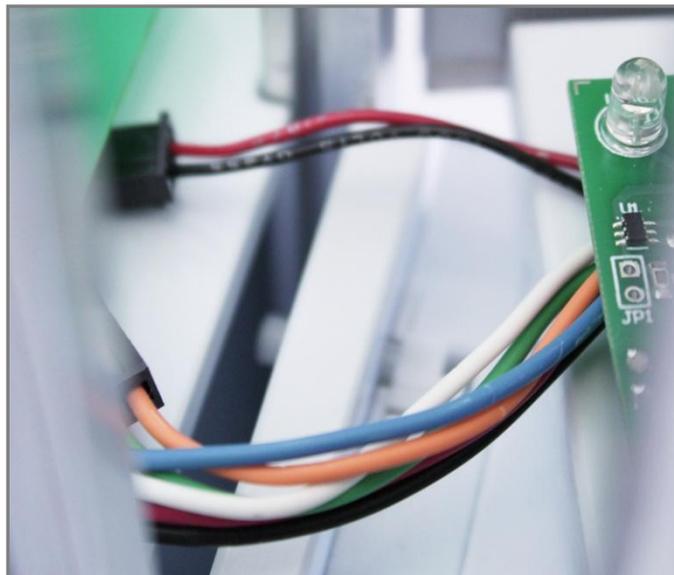

6. Align the snaps on the bottom of the runway with the holes in the hub and press down until they click into place.

*Figure 5 - Runway Snapped into Place*



7. Install the remaining runways in the same manner, connect each cable, tuck the connector and excess cable under the board and snap the runway in place.

## **DISASSEMBLY INSTRUCTIONS**

Firmly grasp the runway and carefully lift up to unsnap it from the hub. Use caution when removing runways; the cables can be pulled if too much force is used.

## ***Appendix A***

### ***Contact Information***

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at [www.med-associates.com](http://www.med-associates.com) for contact information.

For technical questions, email [support@med-associates.com](mailto:support@med-associates.com).