

FEEDER MOUNT FOR MOUSE BACKLIT MAZES

ENV-343U-BLFM

INSTALLATION GUIDE

DOC-257

Rev. 1.0

Copyright ©2014
All Rights Reserved

Med Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478

Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

THIS PAGE INTENTIONALLY LEFT BLANK

Table of Contents

Chapter 1	1
Overview.....	1
Specifications	1
Hardware	1
Chapter 2	2
Assembly.....	2
Mounting the Pellet Feeder	3
Adjusting Feeder Height	4
Appendix A	5
Cleaning.....	5
Appendix B	5
Contact Information	5

CHAPTER 1

Overview

The ENV-343-BLFM feeder mount bracket allows for an optional pellet feeder (ENV-203) to be mounted on a backlit mouse runway. It mounts to the support at the end of the runway.

Specifications


Pellet Feeder Mount Dimensions:

Length:	2.9"
Width:	4.0"
Height:	1.7"

Hardware

The Feeder Mount Bracket holds the Pellet Dispenser at the runway terminus.

Figure 1.1 –ENV-343U-BLFM Feeder Mount with Bracket

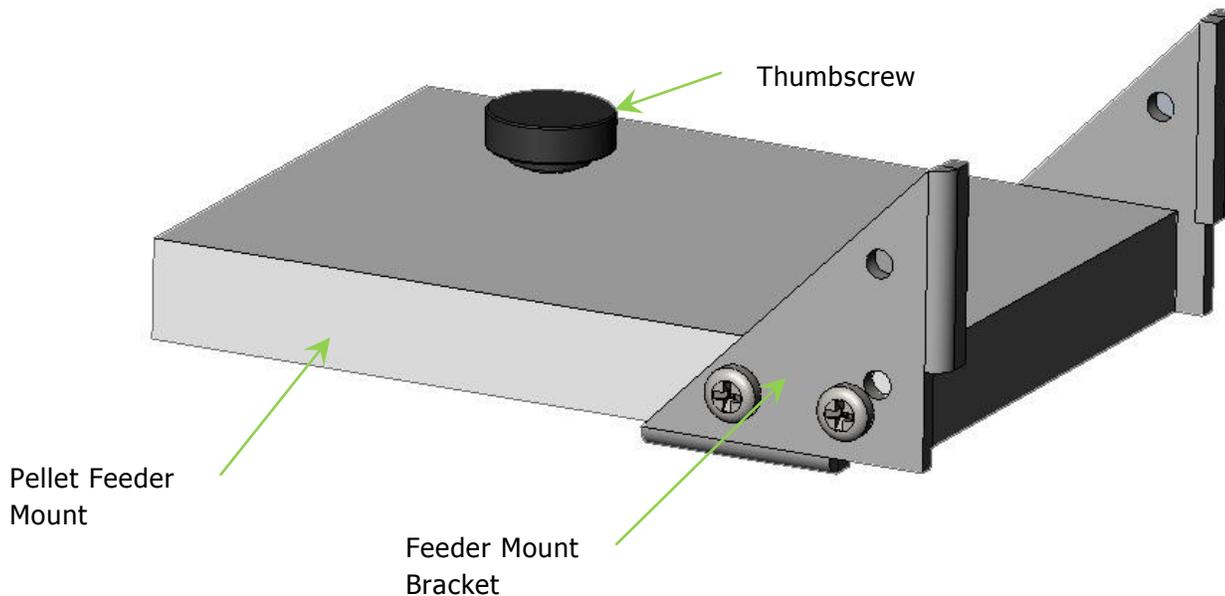

CHAPTER 2

Assembly

Slide the Feeder Mount Bracket onto the runway assembly.

Figure 2.1 –Runway Feeder Mount Bracket Assembly


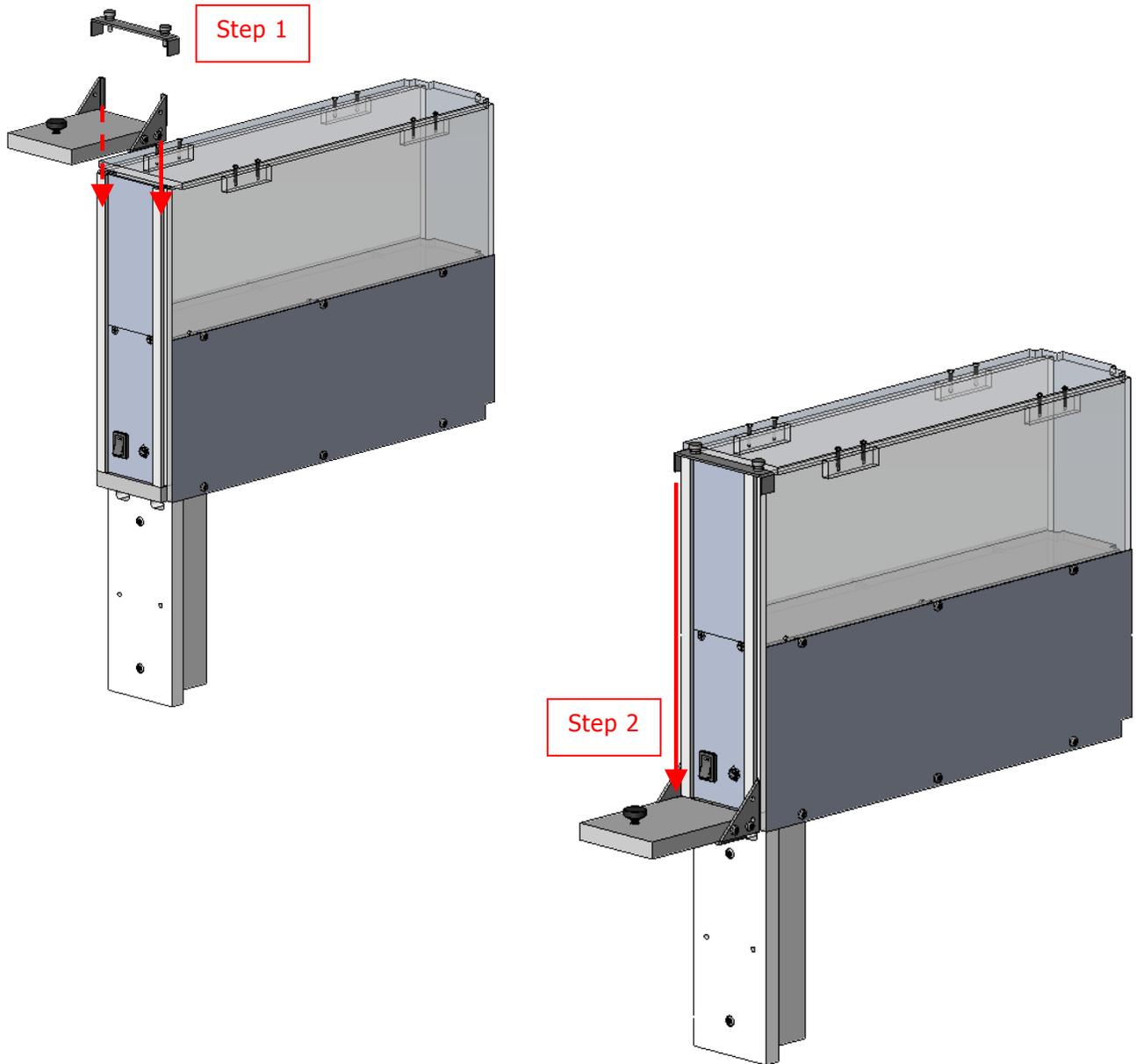


Figure 2.2 – ENV-203 Pellet Dispenser

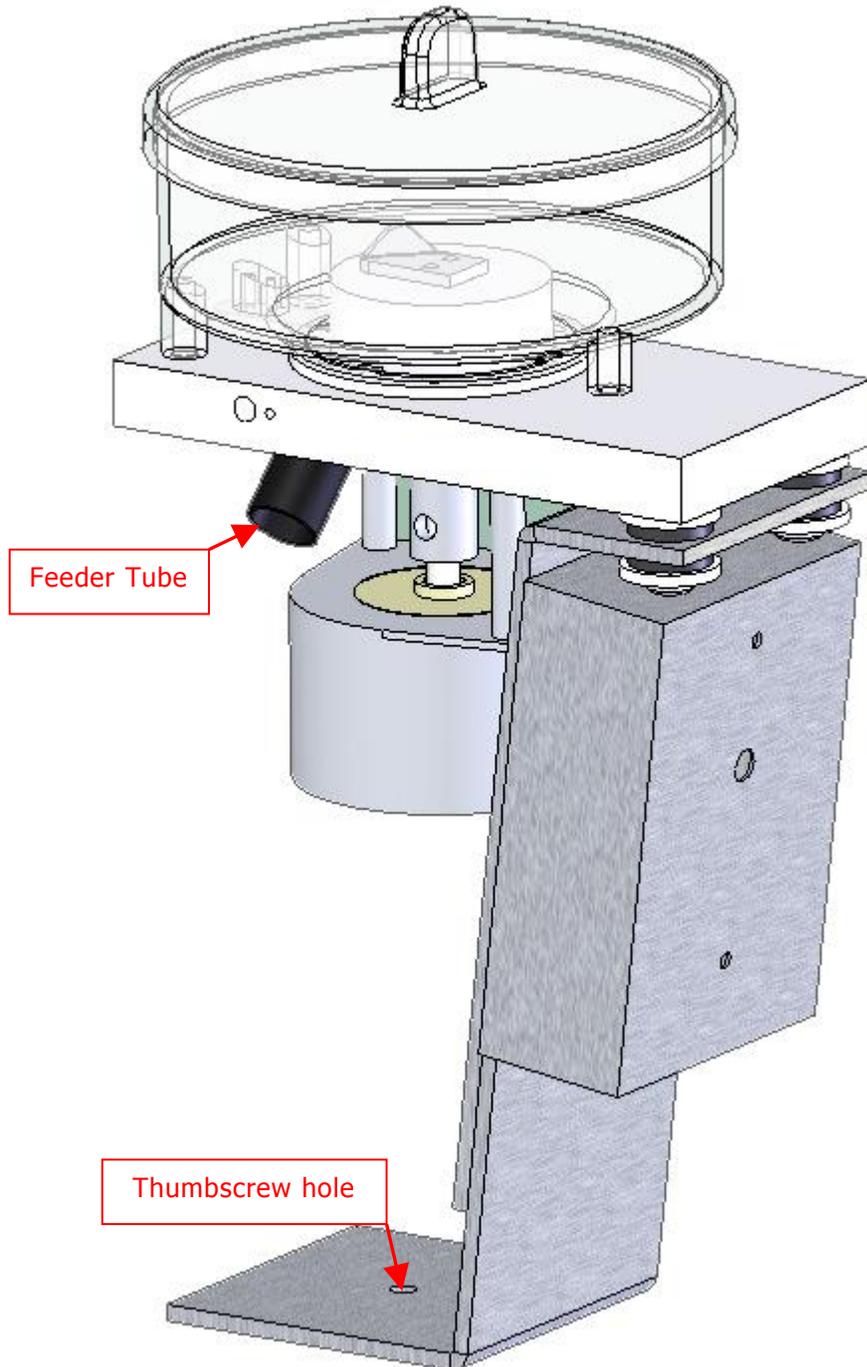


Mounting the Pellet Feeder

Connect the pellet dispenser to the Feeder Mount and secure using the thumbscrew provided. The tube from the pellet dispenser is connected to the tube on the receptacle

with another short piece of tubing (included) so that the pellet is dispensed to the runway.

Figure 2.3 – Final Assembly

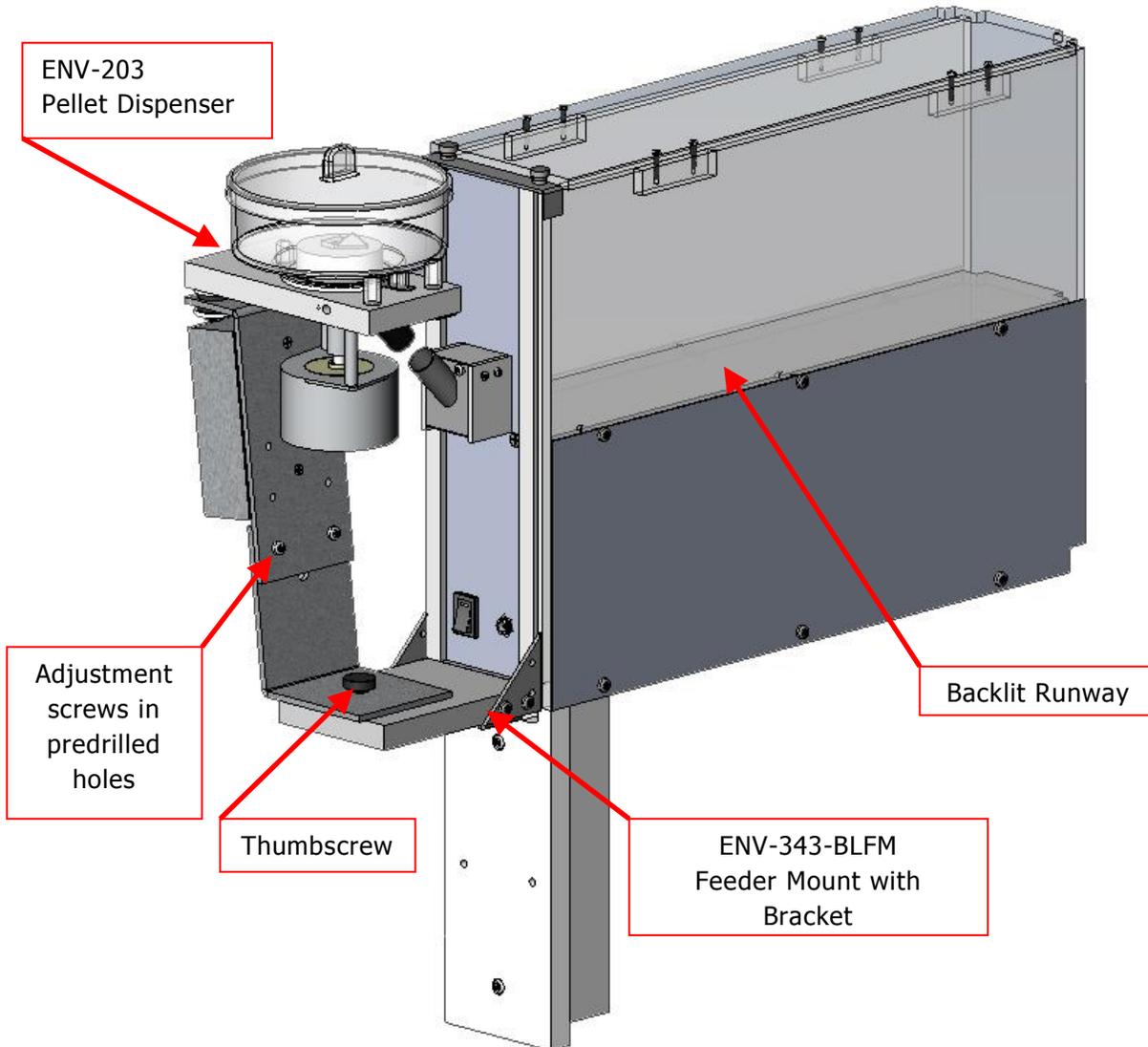

Adjusting Feeder Height

If feeders were purchased separately, the standing height of the feeder may need to be adjusted. If the feeder tube is not aligned with the pellet receptacle chute when the Pellet Dispenser is mounted, remove the adjustment screws and reinstall in the lowest set of holes as shown in Figure 2.3.

Appendix A

Cleaning

Safe for wipe down with Quatricide[®]. A cage washer may be used to clean the feeder mount. Do not exceed 180°.

Appendix B

Contact Information

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.