

A-FRAME INSERT

ENV-008-IRT

USER'S MANUAL

DOC-260

Rev. 1.1

Copyright ©2012
All Rights Reserved

Med Associates Inc.
P.O. Box 319
St. Albans, Vermont 05478

Phone: 802.527.2343
Fax: 802.527.5095
www.med-associates.com

notes

Table of Contents

Chapter 1 | Introduction 1
Specifications 1
Chapter 2 | Hardware 2
Installation 3
Appendix A | Contact Information 4

CHAPTER 1 | INTRODUCTION

The ENV-008-IRT (Rev 1.2) includes a black plastic hinged A-frame insert that is made to fit into chambers ENV-008 and VFC-008. It creates a sloping roof within the chamber from floor to ceiling on two sides and allows unencumbered transmission of NIR light.

Support hardware is included in the ENV-008-IRT to raise the A-frame insert off of the floor. This manual describes how to install the support hardware.

Specifications


Dimensions:

A-Frame Insert	11.6"W x 9.2"D x 7.8"H (29.5 cm x 23.5cm x 19.8 cm)
Support Panel	3.063"W x 3.25"H (7.8 cm x 8.3 cm)

CHAPTER 2 | HARDWARE

The A-frame insert comes with 4 supports, 4 magnets, and 4 support panels. These parts are installed into the chamber to hold the A-frame insert above the floor, allowing room for rats to move under the A-frame insert.

Figure 2.1 – Exploded view of Support Panel and Support

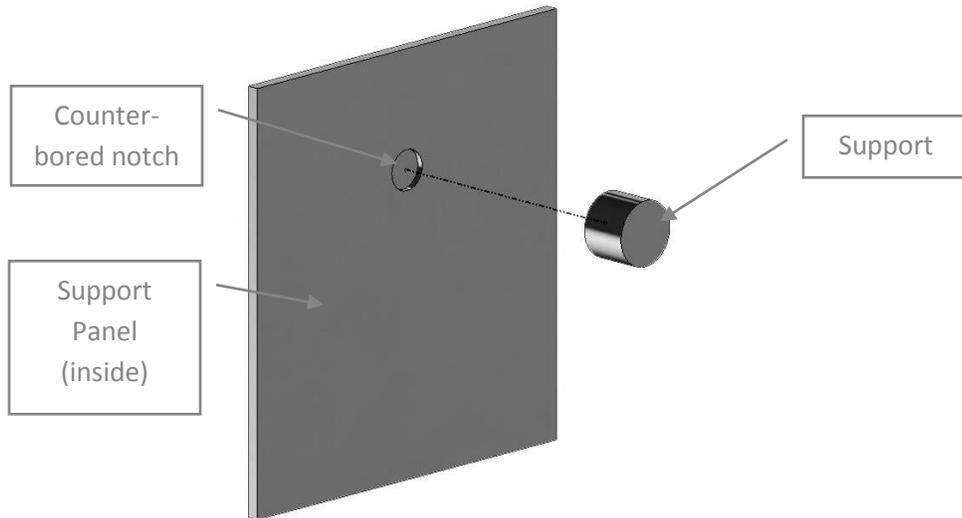

The magnet is placed on the support panel on the outside of the chamber, as shown below.

Figure 2.2 – Magnet Placed on the Support Panel on the Outside of the Chamber


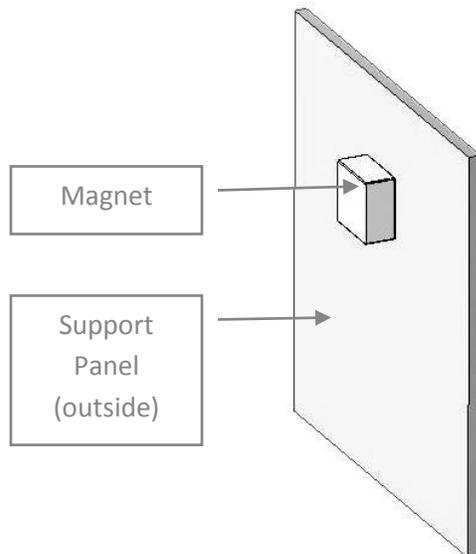


Figure 2.3 –Chambers with A-Frame Insert on Supports

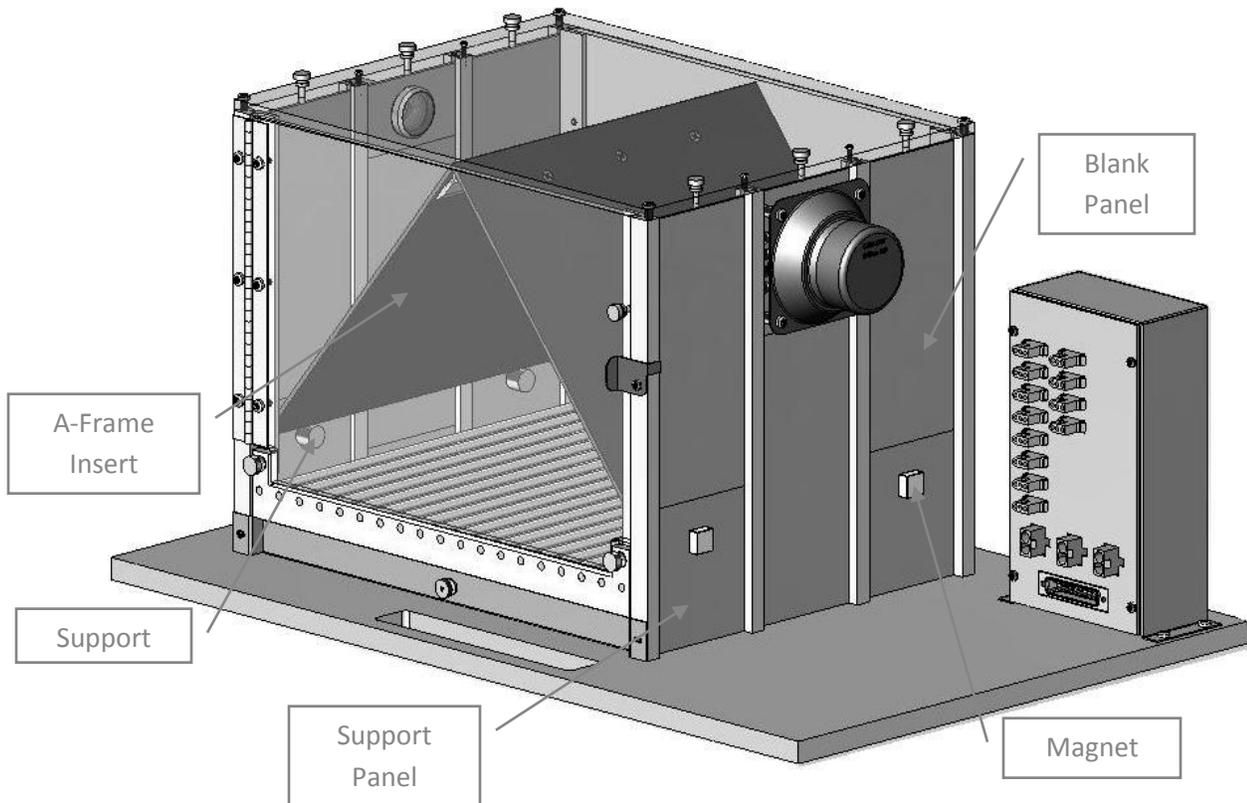

Installation

1. Remove the top of the chamber.
2. Slide out the left and right side blank panels.
3. Insert the support panels.
4. Slide blank panels over the support panels.
5. Place the magnet on the outside of the support panel and the support on the inside of the support panel so that the magnet holds the support. Leave the supports below the counter-bored notch.
6. Repeat for all 4 corners.
7. Place the A-Frame Insert onto the supports.
8. Slide the magnet up so that the support falls into the counter-bored notch in the support panel. The supports are locked into place when you hear a click sound.

Appendix A | Contact Information

Please contact MED Associates, Inc. for information regarding any of our products.

Visit our website at www.med-associates.com for contact information.

For technical questions, email support@med-associates.com.