

# **SOUND ATTENUATING CUBICLE (SAC) REPLACEMENT FAN**

ENV-025F28

ENV-025FAC

INSTALLATION MANUAL

**DOC-313**

**Rev. 1.0**

Copyright ©2017  
All Rights Reserved

Med Associates, Inc.  
P.O. Box 319  
St. Albans, Vermont 05478

Phone: 802.527.2343  
Fax: 802.527.5095  
[www.med-associates.com](http://www.med-associates.com)

## CHAPTER 1 | INSTALLATION

To install the ENV-025F SAC replacement fan, first clip away the back of the white plastic rivet inside the SAC. This will remove the fan from the inside of the SAC and the fan grill from the outside.

Remove screw  
and cable guide.



Align the replacement fan with the cable on the bottom and the side of the fan with the grill inside the SAC. Using the four new plastic rivets secure the grill outside the SAC and the replacement fan inside the SAC. The rivets should press in through the grill on the outside, the SAC wall and the back of the fan on the inside of the SAC when all of the holes are aligned properly. This process may require some force. A rubber mallet may be required.



## APPENDIX A | CONTACT INFORMATION

Please contact Med Associates, Inc. for information regarding any of our products.

For Technical questions, email [support@med-associates.com](mailto:support@med-associates.com).

For Sales questions, email [sales@med-associates.com](mailto:sales@med-associates.com).

Visit our website at [www.med-associates.com](http://www.med-associates.com).